

Features

- Meets Class AAA ASTM standards
- Fiberized output for flexible illumination.
- Easy-to-use touchscreen controls
- Light intensity control is available
- Remote control by PC
- Cost-effective for researchers

Applications

- Testing photovoltaic cell performance
- Photochemistry
- Photobiology
- Testing for color fading of paints and fabrics, etc.





I. Overview


The Sciencetech Solar LightLine series of fiberized solar simulators are a revolution in the solar simulator industry. With the Solar LightLine you can direct the Class AAA solar light anywhere you wish with the standard 4 foot fiber. Classic designs for solar simulators require you to bring your sample to the solar simulator. Now you can bring the solar simulator to your sample.

The Solar LightLine simulators come standard with a touchscreen power supply, integrated shutter, and exposure controller. Multiple filter options are available to meet your specific needs.

The Solar LightLine has four base models (A1-300W, A4-1000W, AX-300W, and CX-300W). Each model has multiple configuration options. Configuration options exist for Class

AAA illumination up to 50mm square and intensity up to 12Suns.

The Sciencetech Solar LightLine can match class A ASTM, IEC, and JIS solar simulator standards. A number of air mass filters are available for the Solar LightLine giving you the capability to customize the instrument for your requirements. In addition to spectral match filters you can also add intensity filters and other spectral filters like UV and IR blocking filters.

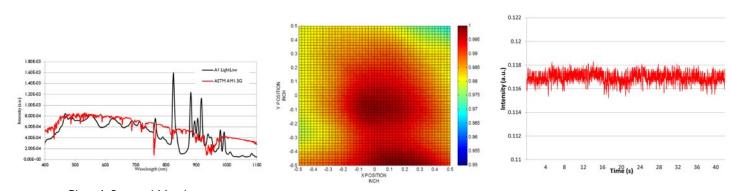

Class A Spectral Match

Class A Non-Uniformity


Class A Temporal Instability

Model	Power operation	Target Size	Sun Level ^l	Working Distance ²	Class ³	Projection Optics
A4-LA125	120/220VAC	25mm x 25mm	Up to 4 Suns	125 to 175 mm	AAA	-LA125
A4-LA200	120/220VAC	50mm x 50mm	I Sun	225 to 300 mm	AAA	-LA200
A4-C250	120/220VAC	Ø12.5mm	Up to 12 Suns	125 to 175 mm	ACA	-C250
AI-LAI25	120/220VAC	25mm × 25mm	I Sun	125 to 175 mm	AAA	-LA125
A1-C250	120/220VAC	Ø12.5mm	4 Suns	125 to 175 mm	ACA	-C250
AX-LA125	120/220VAC	20×20 mm	Up to 3 Suns	125 to 175 mm	AAA	-LA125
AX-LA200	120/220VAC	46×46 mm	I Sun	225 to 300 mm	AAA	-LA200
CX-C250	120/220VAC	Ø12.5mm	9 Suns	125 to 175 mm	ACA	-C250

- I With AMI.5G filter installed
- 2 Working distance can be customized with different optics but will change target size
- 3 Tested to ASTM E927-05 standards, other standard testing is available upon request



2. Specifications Al LightLine

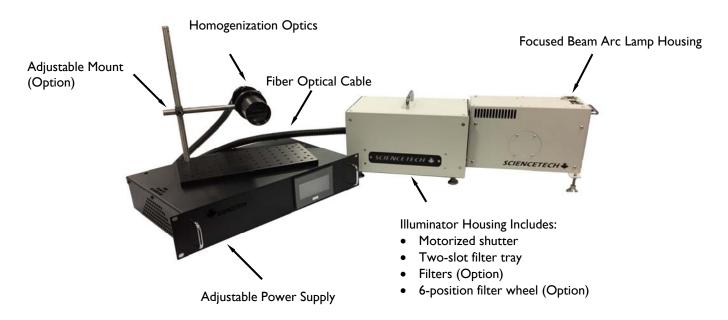

MODEL	AI-LAI25	A1-C250	
ASTM Classification	AAA	ACA	
Target Size	25mm × 25mm	Ø12.5mm	
Sun Level	l Sun	4 Suns	
Working Distance	125 to 175 mm	125 to 175 mm	
Wavelength Range	350 - 1800nm (Standard Spectral Match 400 - 1100nm)		
Air Mass Filter	AM1.0D, AM1.5G, AM1.5D, AM2.0		
Collimation Angle	10° Half Angle (depends on selected optics)		
Lamp	300W Xenon, 1000h (average) Lifetime, Forced Air Cooling		
Shutter	Integrated Touchscreen Control, 0 - 10000s, Adjustable Duty Cycle		
Fiber Bundle Length	I.22m (4') Standard		
Power Supply Model	601-300		
Input Voltage/Current	Switchable I I0-I 30V / 3A or 220-240V / I.5A Operation		
Environment	Temperature 10-35 °C, Humidity 20-80%, Avoid Condensation		
Dimension (LxWxH)	58.7cm × 20.6cm × 20.8cm		
Weight	5kg		



LightLine X




MODEL	AX-LA125	AX-LA200	CX-C250
ASTM Classification	AAA	AAA	ACA
Target Size	20x20 mm¹	46×46 mm ¹	Ø12.5mm
Sun Level	3 Suns	I Sun	9 Suns
Working Distance	125-175 mm	225-300 mm	125-175 mm
Wavelength Range	350 - 1800nm (Standard Spectral Match 400 - 1100nm)		
Air Mass Filter	AMI.0D, AMI.5G, AMI.5D, AM2.0		
Collimation Angle	10° Half Angle (depends on selected optics)		
Lamp	300W Xenon, 1000h (average) Lifetime, Forced Air Cooling		
Shutter	Integrated Touchscreen Control, 0 - 10000s, Adjustable Duty Cycle		
Fiber Bundle Length	I.22m (4') Standard		
Power Supply Model	601-300		
Input Voltage/Current	110-130V / 3A or 220-240V / 1.5A		
Environment	Temperature 10-35 $^{\circ}\text{C}$, Humidity 20-80% , Avoid Condensation		
Dimension (LxWxH)	76.8cm × 21.6cm × 38.3cm		
Weight	20kg		

 $^{^{\}rm I}$ Up to 25% larger classified uniform areas available at Class B and C ASTM non-uniformity.



A4 LightLine

Illuminator Housing Included:

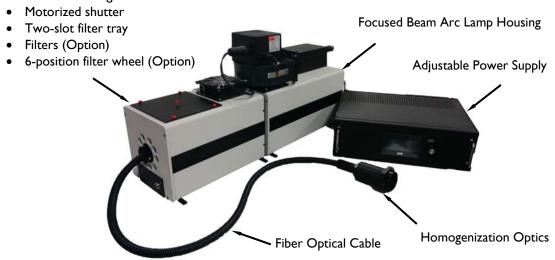

MODEL	A4-LA125	A4-LA200	A4-C250
AST Classification	AAA	AAA	ACA
Target Size	25mm × 25mm	50mm x 50mm	Ø12.5mm
Sun Level	Up to 4 Suns	I Sun	Up to 12 Suns
Working Distance	125 to 175 mm	225 to 300 mm	125 to 175 mm
Wavelength Range	350 - 1800nm (Standard Spectral Match 400 - 1100nm)		
Air Mass Filter	AMI.0D, AMI.5G, AMI.5D, AM2.0		
Collimation Angle	10° Half Angle (depends on selected optics)		
Lamp	1000W Xenon, 1000h (average) Lifetime, Forced Air Cooling		
Shutter	Integrated Touchscreen Control, 0 - 10000s, Adjustable Duty Cycle		
Fiber Bundle Length	I.22m (4') Standard		
Power Supply Model	6II-Ik		
Input Voltage/Current	110-130V / 9A or 220-240V / 6A		
Environment	Temperature 10-35 °C , Humidity 20-80% , Avoid Condensation		
Dimension (LxWxH)	76.8cm × 21.6cm × 38.3cm		
Weight	20kg		




3. Accessories

Touchscreen Power Supply

MODEL (Default)	DESCRIPTION
Shutter and Exposure Controller	This feature provides an onboard shutter and exposure controller built into the power supply. A standard 2-pin connector is provided on the back of the power supply to connect a Sciencetech solenoid driven shutter mechanism. With this option the power supply firmware provides an interface for programming shutter loops and exposure sequences.
Power Connection for Cooling Fans	The power connection for cooling fans feature provides a power pass through connection for the lamp house cooling fans. With this option the lamp housing cooling fans operation is monitored by the lamp house.
Safety Interlock	With the safety interlock feature the power supply will turn off, extinguishing the lamp, if the interlock is opened. The interlock wiring can be connected to a mechanical safety interlock such as a magnetic switch or rocker switch on the lamp house.
RS232 Computer Control	This feature provides RS232 computer control of the power supply through a DB9 connector mounted on the rear panel of the power supply (Power control 1.1 software and cable are included).
MODEL (Option)	DESCRIPTION
Programmable Remote Temperature Monitor/TM	The programmable remote temperature monitor option provides over temperature protection for the lamp house that the power supply is being used with. With this option an RTD or K-type Thermocouple is wired between the lamp house and power supply.
Monitor Output/OC	This option provides a male nine pin connector on the rear of the power supply for monitoring the status of the power supply. Parameters that can be monitored: lamp current 0-10V, lamp voltage 0-10V, lamp On/Off (1-5V TTL).
Auto Start Lamp/AS	The option provides the capability to automatically start the lamp. This allows the user to set a time for the light source to automatically ignite.



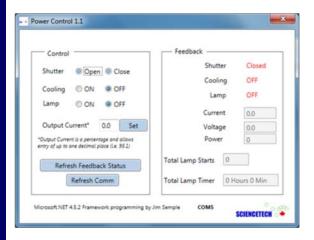







Filters

Numerous filtering options are available with the Sciencetech Solar LightLine. The standard Solar LightLine comes with two filter slots. One slot is usually reserved for an air mass filter. The secondary slot can accommodate other filtering options.

MODEL	Code	DESCRIPTION
FT-75-3	100-8035	Empty FT Style Filter Holder for 75mm Square or Diameter Filters (3mm thick)
MF-34-FT-3	640-9005	75mm Mesh 34% Neutral Density Filter with Holder
MF-49-FT-3	640-9006	75mm Mesh 49% Neutral Density Filter with Holder
MF-71-FT-3	640-9007	75mm Mesh 71% Neutral Density Filter with Holder
HPF-BB-NIR-FT-3	640-9012	Passes UVA, UVB, and VIS (280-730nm). Blocks NIR (730-1100nm).
HPF-LP-VIS-FT-3	640-9014	Passes VIS and NIR. Blocks UVA and UVB (<400nm).
HPF-LP-NIR-FT-3	640-9015	Passes NIR (700-2000 nm). Blocks UVA, UVB, and VIS (<700 nm).
HPF-BB-VIS-FT-3	640-9016	Passes UVA and UVB (245-400nm). Blocks VIS (400-700nm).
HPF-LP-UVA-FT-3	640-9017	Passes UVA,VIS, and NIR. Blocks UVB (<320nm)
HPF-BB-UVB-FT-3	640-9018	Passes UVB, VIS, and NIR. Blocks UVA (320-400nm)
HPF-BP-VIS-FT-3	640-9019	Passes VIS (400-700nm). Blocks UVA, UVB, and NIR.
MF-25	640-9013	Set of 25mm Diameter Wire Mesh Filters to Attenuate Solar Level. For use with the –SFW option.
AMI.0D-FT-3	160-8022	Air Mass AMI.0D (Standard Range)
AMI.5G-FT-3	160-8024	Air Mass AM1.5G (Standard Range)
AMI.5D-FT-3	160-8026	Air Mass AM1.5D (Standard Range)
AM2.0D-FT-3	160-8028	Air Mass AM2.0D (Standard Range)
WF-1Q	100-8048	Compact IR Water Filter



Mesh filter











Optional Accessories

MODEL	PART No.	DESCRIPTION
SC-12	130-9011	A general-purpose sample chamber that can be configured in multiple ways to address a number of application requirements. The SC-12 includes 4 optical ports (on three sides and one on the top) as well as a side door for easy access to the sample area. Three of the side walls feature optical ports with a Sciencetech standard B4 hole pattern as well as removable anodized access panels that can be machined to accommodate most types of feedthroughs, including electrical, liquid and gas.
-PM90	100-8060	A stand/fiber holder designed to hold the LightLine optical illumination delivery fiber. Fiber holder position is adjustable for vertical and horizontal illumination and any angle in-between. Includes a sturdy baseplate.
FB-0.5-8	670-0103	8 foot optical fiber upgrade to allow for further versatility with the A1, AX/CX and A4 LightLine. Please note that there will be a reduction in the power output with an 8 foot fiber.
-SFW	189-9012	6-position computer controlled filter wheel (25 mm diameter filters)
(UV-Glasses-Drk)	720-0159	Dark safety glasses






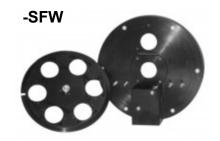


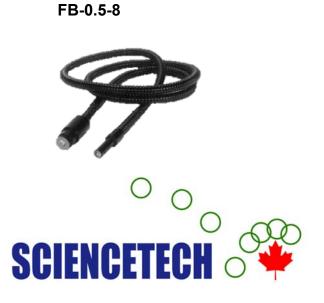


5. Dimensions

AI/AX/CX LightLine

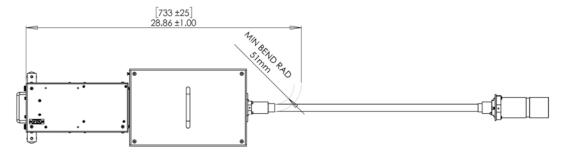

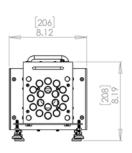

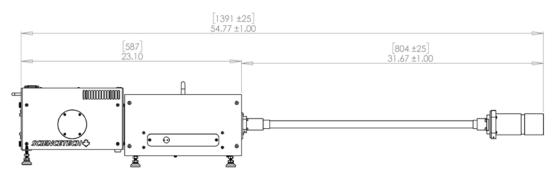

A4 LightLine

